



Sterile ready to use solution based on non-ionic surfactants for the cleaning of surfaces in cleanrooms.

perform® sterile cleaner ND

Our Plus

- ready to use
- sterile (X-Ray irradiated) and microbial filtrated (0.2 µm)
- available in convenient sterile 500 ml spray system
- colour and perfume free
- excellent cleaning performance

Application areas

Cleaning of surfaces and equipment / devices in cleanrooms.

Especially for Grade A and B clean areas, according to:

- EU Guidelines to Good Manufacturing Practice for medicinal products, Annex 1: Manufacture of sterile medicinal products, March 2009.
- FDA Guidance for Industry: Sterile drug products produced by aseptic processing - current Good Manufacturing Practice, U.S. Department of Health and Human Services; Food and Drug Administration, September 2004.

Instructions for use

Apply perform® sterile cleaner ND undiluted to objects and surfaces and wipe if necessary.



perform[®] sterile cleaner ND

Product data

Labelling according to Regulation (EC) No. 648/2004: < 5 % non-ionic surfactants.

Chemical-physical data

Color	colourless
Density	approx. 1,00 g/cm ³ / 20 °C
Flash point	Not applicable
Form	liquid
pH	approx. 5,8 / 20 °C
Viscosity, dynamic	not determined

Special advice

The sterile 500 ml spray system offers the following advantages:

- convenient and space saving
- ergonomically shaped spray pump will not slip out of a gloved hand
- still sterile after opening

The bottles are sterile (with radiation indicator) and double-wrapped individually.

Information for order

Item	Delivery form	Item no.
perform [®] sterile cleaner ND 500 ml bottle	10/Carton	on request

These products are not available in every country. For more information please contact our local subsidiary or distributor.

Related Products

- perform[®] sterile dry wipes
- perform[®] sterile mix dry wipes
- perform[®] sterile premium dry wipes

Environmental information

schülke manufactures products economically and with advanced, safe and environmentally friendly production processes while at the same time maintaining our high quality standards.

Expert opinion and information

Please visit our website for an overview of all available literature/reports on the product: www.schuelke.com.

For individual questions:

Customer Sales Service

Phone: +49 40 52100-666

E-Mail: info@schuelke.com



Schülke & Mayr GmbH holds a Manufacturer's Authorisation according to sect 13 para 1 German Drug Law and Certificates of GMP Compliance for medicinal products.

schülke Headquarters
Schülke & Mayr GmbH
Robert-Koch-Str. 2
22851 Norderstedt, Germany
Phone +49 40 - 52100 - 0
Fax +49 40 - 52100 - 318
www.schuelke.com
info@schuelke.com